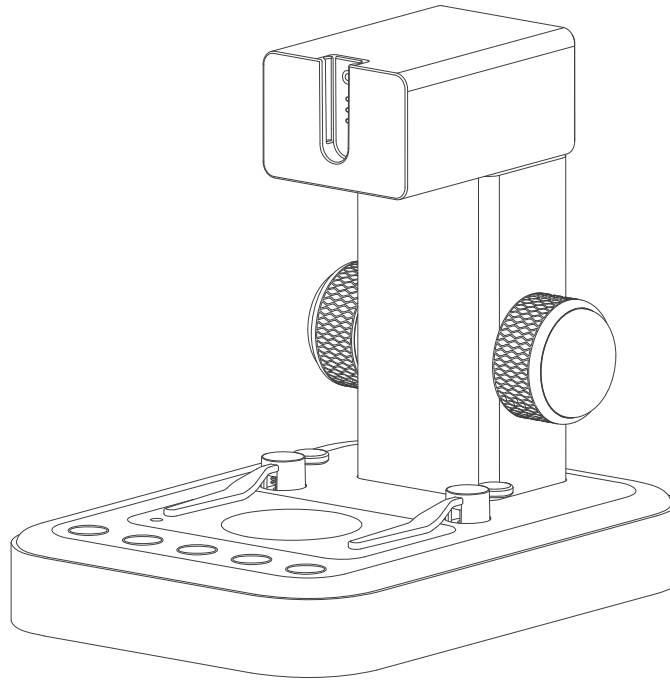
For Digital Microscope MX Series

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MX-Stand

Product description

Thank you for purchasing the MatataStudio MX series stand. This stand features built-in lighting and supports intelligent staining functionality. When used with a microscope, it allows for multi-level hybrid focusing and stepless zooming. The anti-shake design ensures more stable imaging.

Contents EN

Microscope and MX-Stand Installation
 Instructions 1
 MX-Stand Usage Instructions 2
 Steps to Use 3
 Specifications 4
 Parts list 4
 Warranty terms 4

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

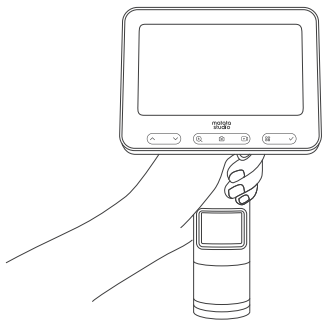
Microscope and MX-Stand Installation Instructions

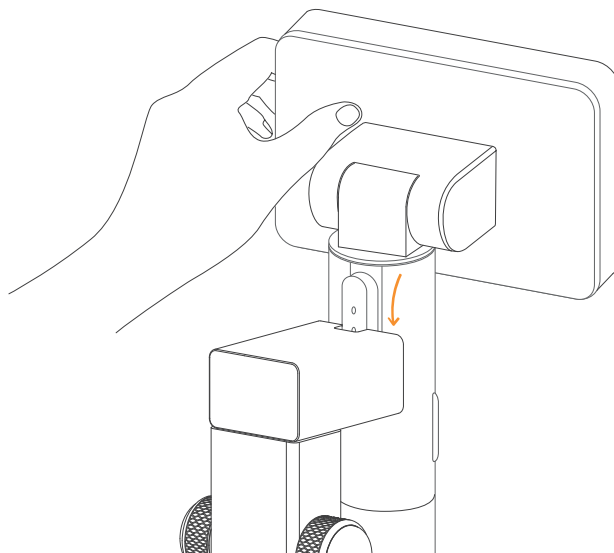1. Hold the microscope steady with one hand.

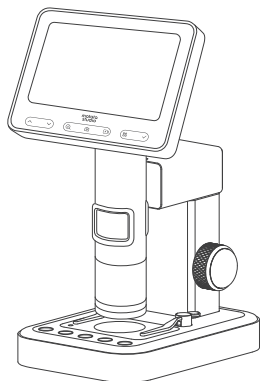
2. Hold the stand with another hand and insert the microscope vertically downward into the mounting hole, making sure the slot at the rear are aligned with the mounting hole.

3. Insert it into the base.

Note: The Stand itself does not contain a power supply. Once the MX-Stand is installed, the stand will be powered when the microscope is turned on.

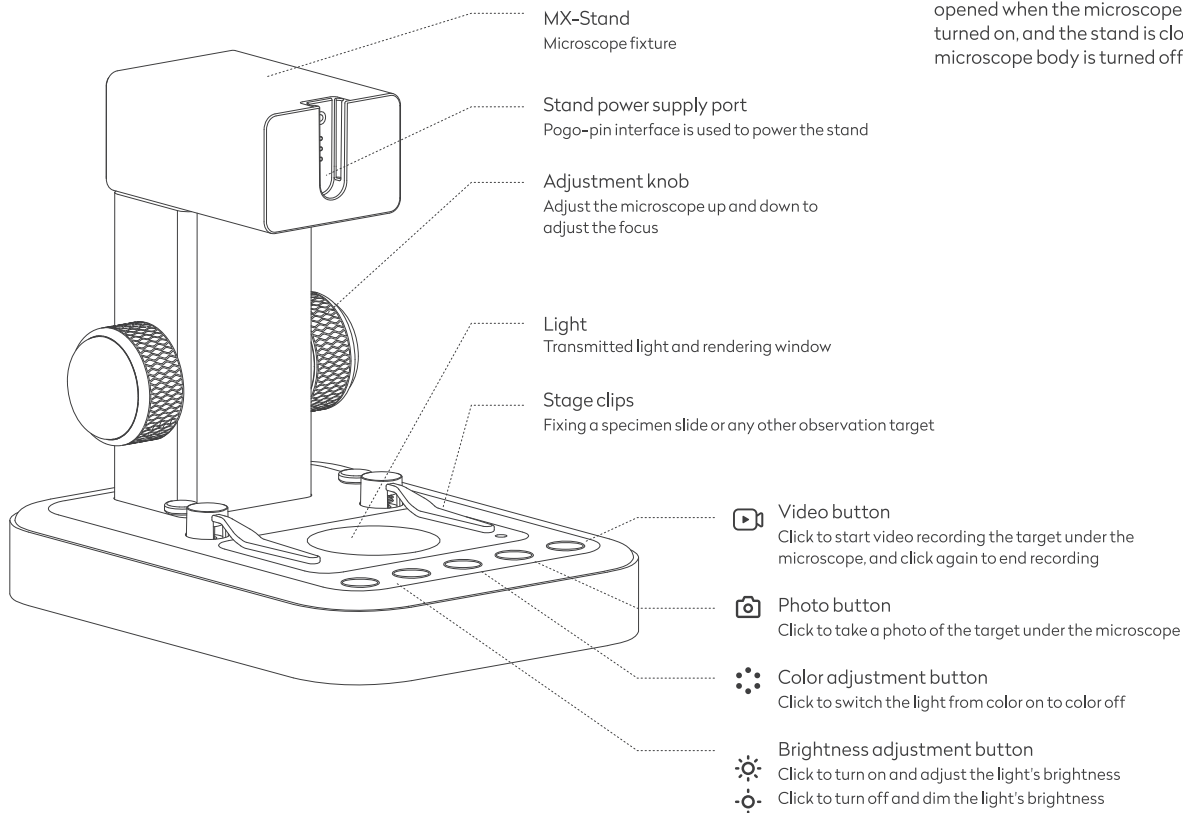
MX-Stand Usage Instructions


Turn on/ off MX-Stand

After the MX-Stand is installed, the stand is turned on when the microscope body is turned on, and the stand is closed when the microscope body is turned off.

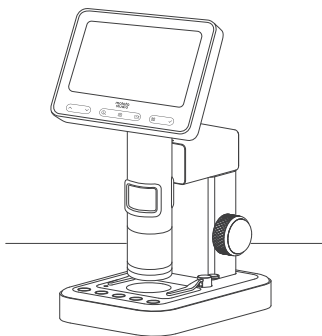
Steps to Use

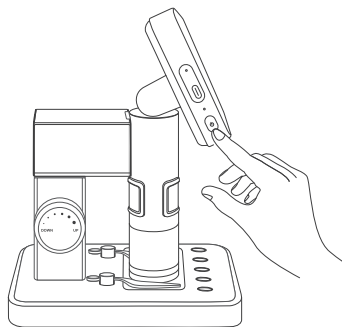
1. Place the microscope on a stable table to avoid shaking or falling.

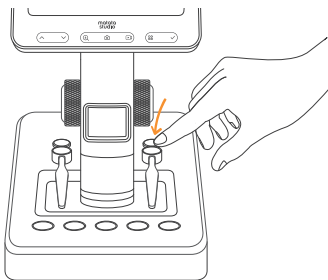
2. Turn on the microscope body.

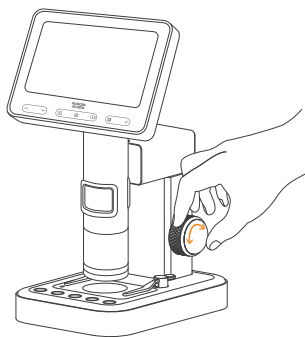3. Place the object to be observed in the center of the bottom stage.

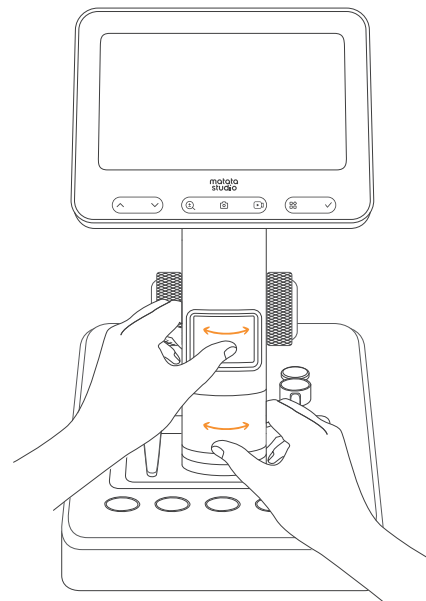
4. Adjust the microscope lifting knob to make the tube at a relatively appropriate height for easy observation of the object.


5. Adjust the focus and objective lens knob, observe the screen display, and slowly adjust for the best observation effect.

Note: You can adjust the observation height and focal length to achieve different magnifications of observation.

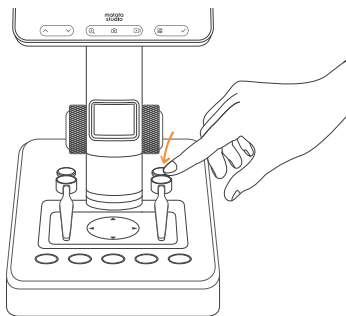Use of stage clips

Use of the stage

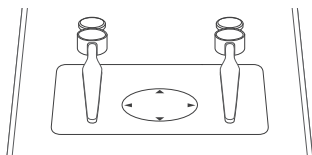Note: Please wipe the stage regularly to avoid affecting the observation experience.


Description of the light adjustment function

- The intensity of the light will affect the specimen's outline and some details. Please pay attention and adjust as needed during use.
- Choose the right rendering light to achieve the best observation effect of the specimen.


Specifications

| | |
|---------------------------|-------------------------|
| Length, width, and height | 150.5mm*120mm*168mm |
| Material | Aluminum alloy exterior |
| Weight | 480g |

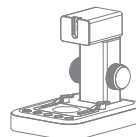Parts list

MX-Stand

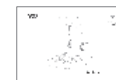User guide

Warranty terms

*Product warranty period is 1 year

*The following situations are not covered by free warranty:

- Unable to present this warranty certificate and valid invoice
- This warranty certificate contains traces of alteration or is inconsistent with the product
- Natural consumption/wear and aging of consumable materials
- Damage caused by lightning strikes or other electrical system reasons
- Damage caused by improper use by users, such as liquid injection/external force damage, etc.
- Damage caused by force majeure factors such as accidents/disasters
- Products that have been disassembled/modified/repared by yourself
- Damage caused by exceeding the warranty period or not using/maintaining/storing the product in accordance with the product instructions

This device complies with the essential requirements and other relevant provisions of the Low Voltage Directive 2014/35/EU, the EMC Directive 2014/30/EU, the Eco-Design Directive 2009/125/EC and the ROHS Directive 2011/65/EU.

WASTE ELECTRICAL AND ELECTRONIC EQUIPMENT (WEEE)

The WEEE marking indicates that this product should not be disposed with regular household waste at the end of its life cycle. This regulation is created to prevent any possible harm to the environment or human health. This product is developed and manufactured with high quality materials and components which can be recycled and/or reused. Please dispose this product at your local collection point or recycling centre for electrical and electronic waste. This will make sure that it will be recycled on an environmentally friendly manner, and will help to protect the environment in which we all live.

This user guide contains important information, please keep it!

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MATATALAB CO., LTD. in China